

Abaxis Fuse

30/07/2025 11:49 am BST


Summary

The Abaxis Vet Scan Fuse integration allows the ordering of tests from within Merlin directly to the Abaxis lab equipment, when results are available they are automatically saved to the patients history.


Details

How do I create a Lab Request?


1. Select a Patient and enter the Lab Screen.

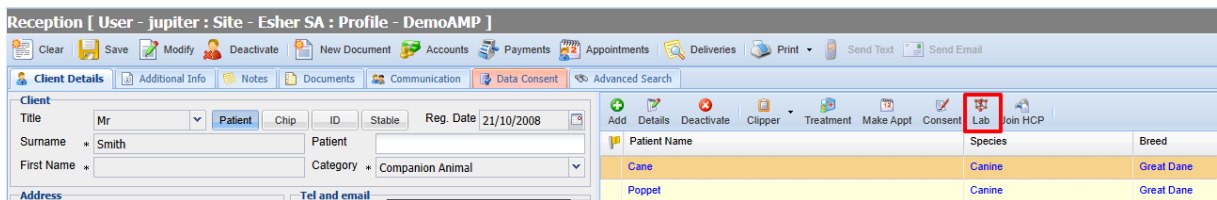


2. Select the 'Vet Scan Fuse' tab.

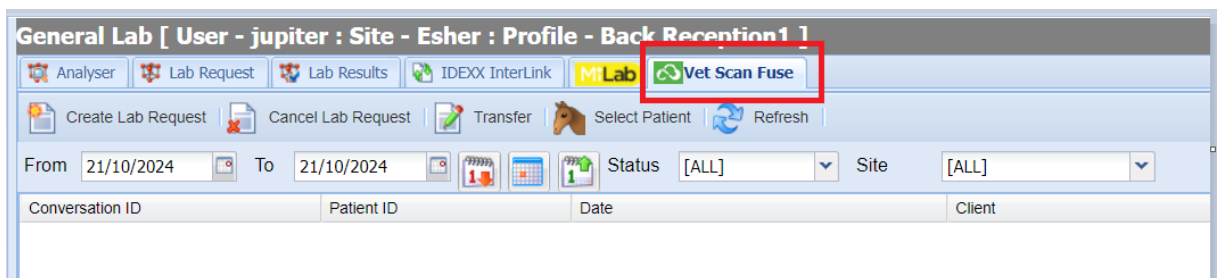

3. Then, select 'Create Lab Request'.

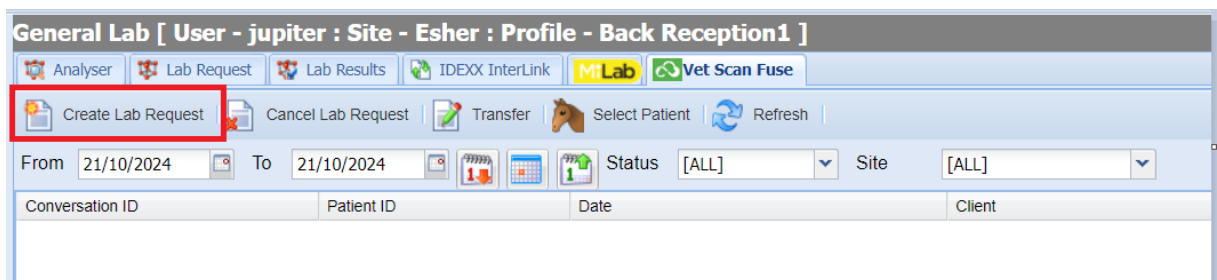

4. The following window will appear. From this window, tick the required lab requests.



Vet Scan Fuse New Order

Create Request Exit Note: Hematology machines cannot accept other species

Sites: Esher

Surgeon: All

Vet Scan Gender:

Vet Scan Species:

Breed: Angora

Lab Requests


Chemistry

☐ Avian/Reptilian Profile Plus(AR)

☐ Canine Wellness Profile(CWP)

☐ Comprehensive Diagnostic(CDP)

5. When selecting a request, you are prompted to select the relevant Abaxis Device.

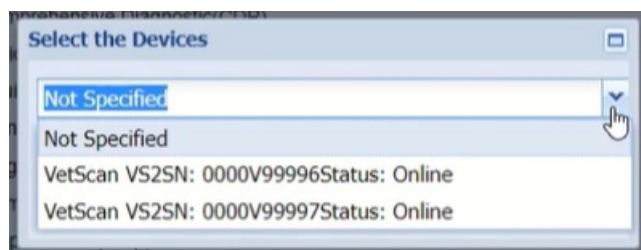

Select the Devices


Not Specified

Not Specified

VetScan VS2SN: 0000V99996Status: Online

VetScan VS2SN: 0000V99997Status: Online


6. Once you have selected the necessary requests and selected the relevant devices, select 'Create Request' to generate.

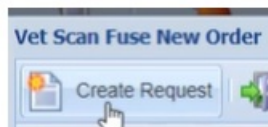

Vet Scan Fuse New Order

Create Request


7. The lab request will display under the Vet Scan Fuse Tab. It will remain under this tab with the updated status displayed in the 'Status' column.

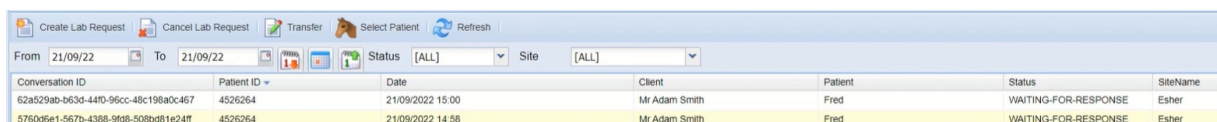

Conversation ID	Patient ID	Date	Client	Patient	Status	SiteName
62a529ab-b63d-44f0-96cc-48c198a0c467	4526264	21/09/2022 15:00	Mr Adam Smith	Fred	WAITING-FOR-RESPONSE	Esher
5760d6e1-567b-4388-9f08-508bd81e24ff	4526264	21/09/2022 14:58	Mr Adam Smith	Fred	WAITING-FOR-RESPONSE	Esher

8. You can double click on the request to view details.

9. From this screen you can cancel the request or print the lab request.

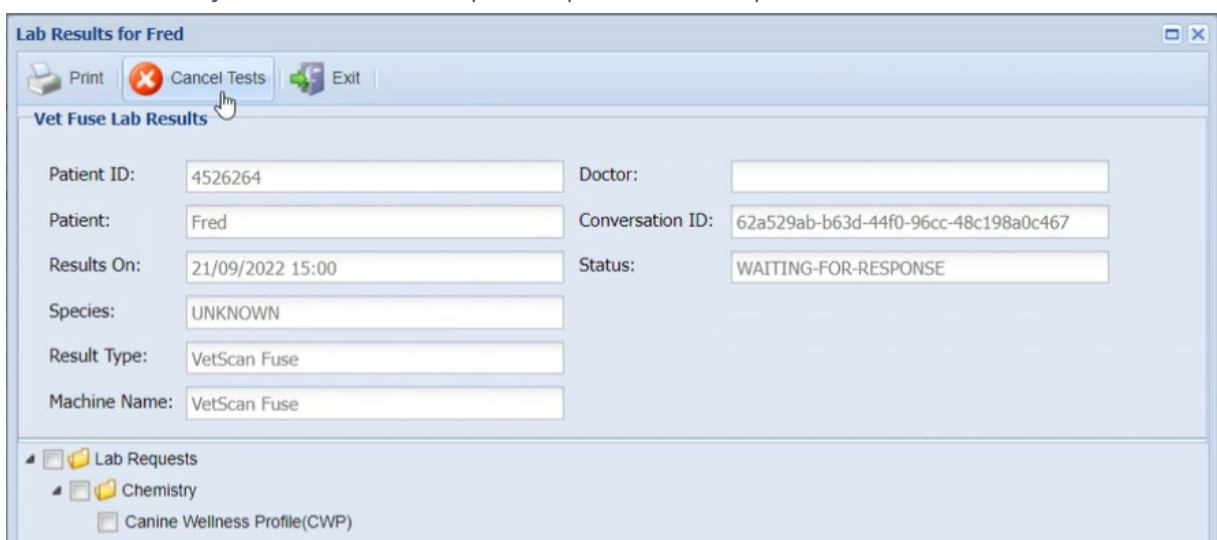

Lab Results for Fred

Print Cancel Tests Exit

Vet Fuse Lab Results

Patient ID: 4526264 Doctor:

Patient: Fred Conversation ID: 62a529ab-b63d-44f0-96cc-48c198a0c467

Results On: 21/09/2022 15:00 Status: WAITING-FOR-RESPONSE

Species: UNKNOWN

Result Type: VetScan Fuse

Machine Name: VetScan Fuse

Lab Requests

Chemistry

☐ Canine Wellness Profile(CWP)

NOTE: any lab requests done directly from the VetScan machine will still show in the results list but they will not be saved in the patients history. To do this click on the test result, select the Transfer button, search for the relevant client and patient and press transfer.



How do I view Lab Results?

Results will be displayed under the Vet Scan Fuse Tab. These can be accessed by double clicking the relevant record.

Alternatively, the results can be accessed via the treatment history screen. When a lab request is created this will display in the treatment screen and when results are available these will display and can be opened by double-clicking the record.

How to Setup Integration

Please contact your Regional Customer Manager (RCM) and our installations team will assist with the configuration with your equipment.
